

# BIOREACTORS

## NEW SOLUTIONS FOR OLD PROBLEMS

Dr. Pavel Lehky

LAMBDA – Laboratory Instruments

[www.rhone.ch/lambda](http://www.rhone.ch/lambda)


# OUTLINE

- Stirring mechanism
- IR-heating system
- Approach used in the MINI FOR
- Probes
- Fermentation accessories (Pumps and gas flow controller)
- Conclusions

# TRADITIONAL STIRRING

## a) Mechanical seals


- No physical separation
- Contaminations
- Service "sensitive"

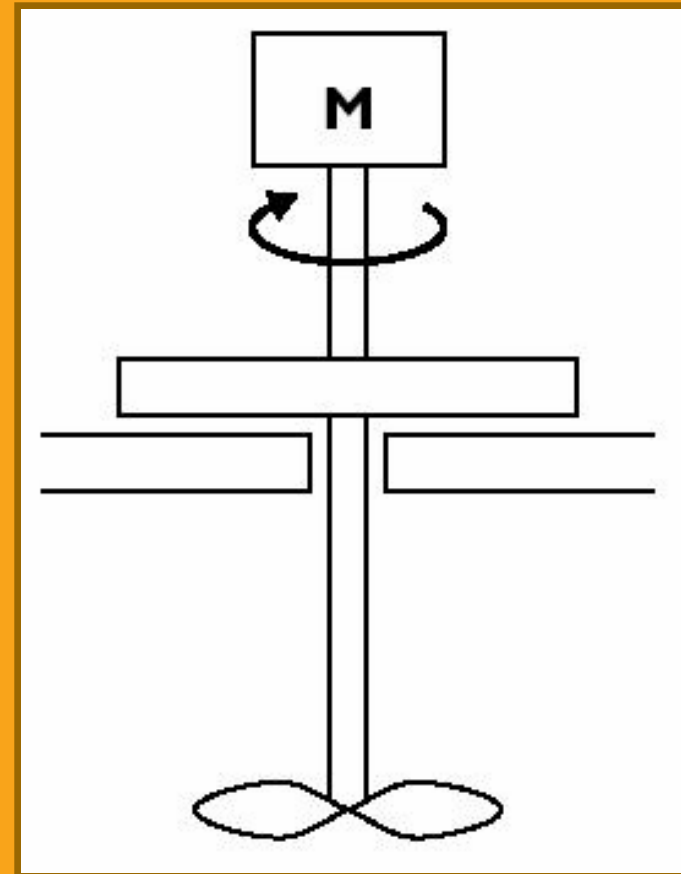

# TRADITIONAL STIRRING


## b) Magnetic coupling

- Separation  $\Rightarrow$  Sterility
- Dead spaces
- High cost

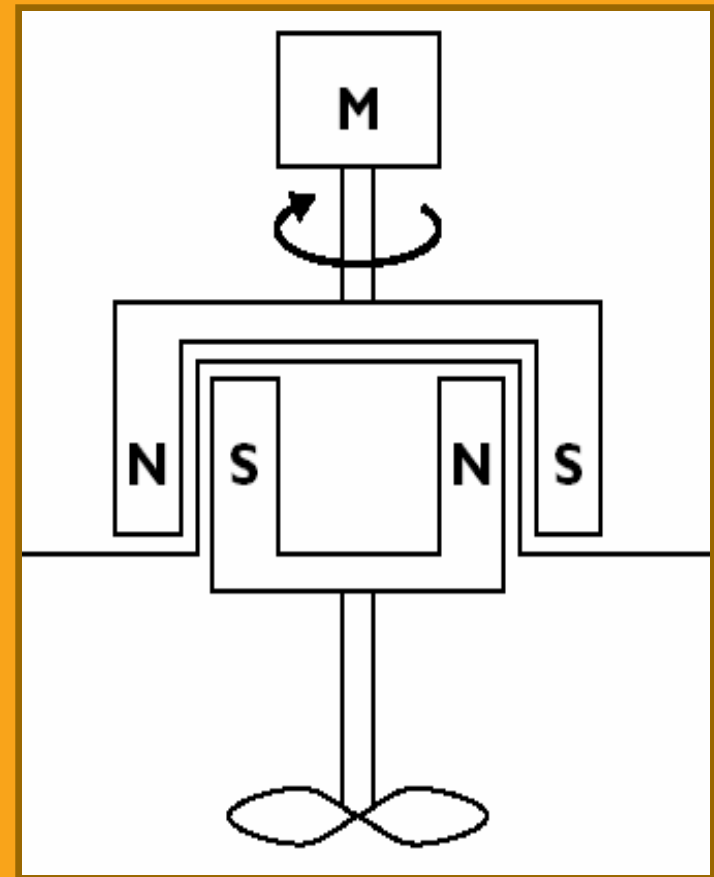

# NEW STIRRING SCHEME


Up and down movement

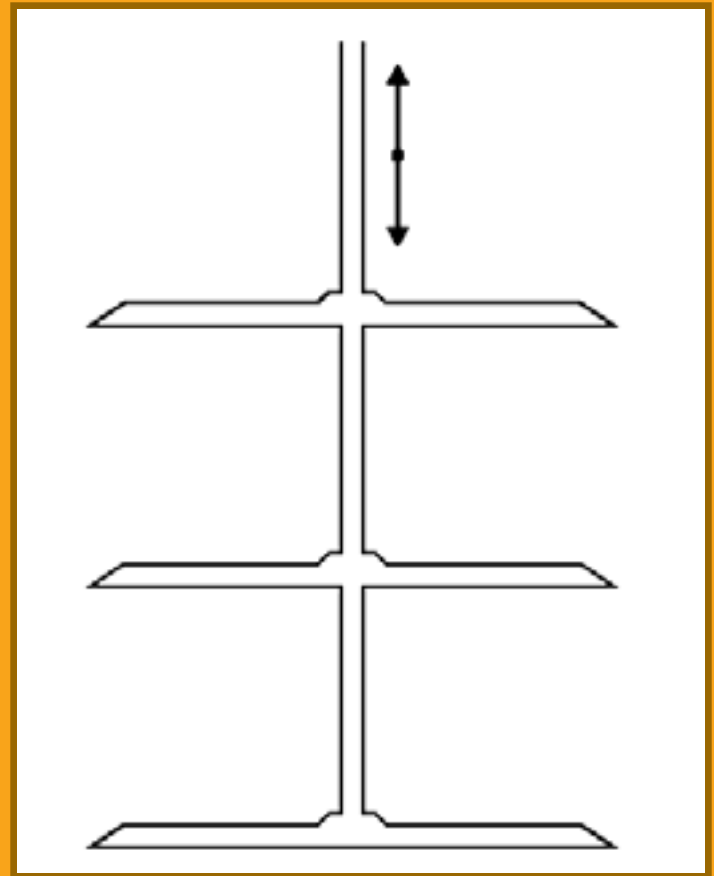

# NEW STIRRING SCHEME

## Fermentor head



- Separation  $\Rightarrow$  Sterility
- Electromagnet  $\Leftrightarrow$  Motor
- Easy to mount (one hand)
- Threaded necks
- Low cost

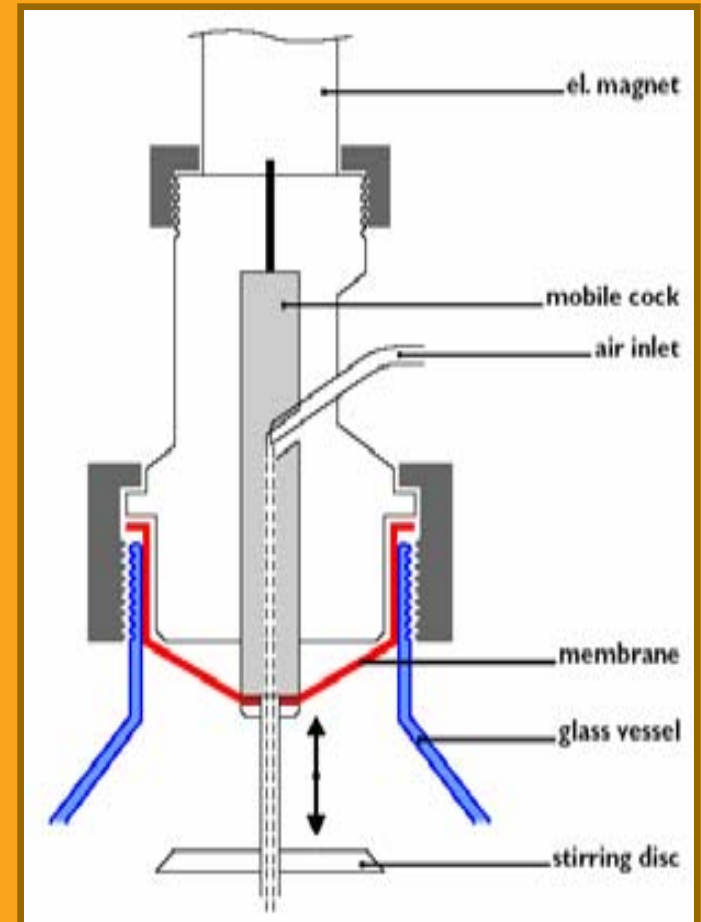

# NEW STIRRING SCHEME

## Stirring discs

- Counter rotating flow paths
- Vertical stream (through the holes)

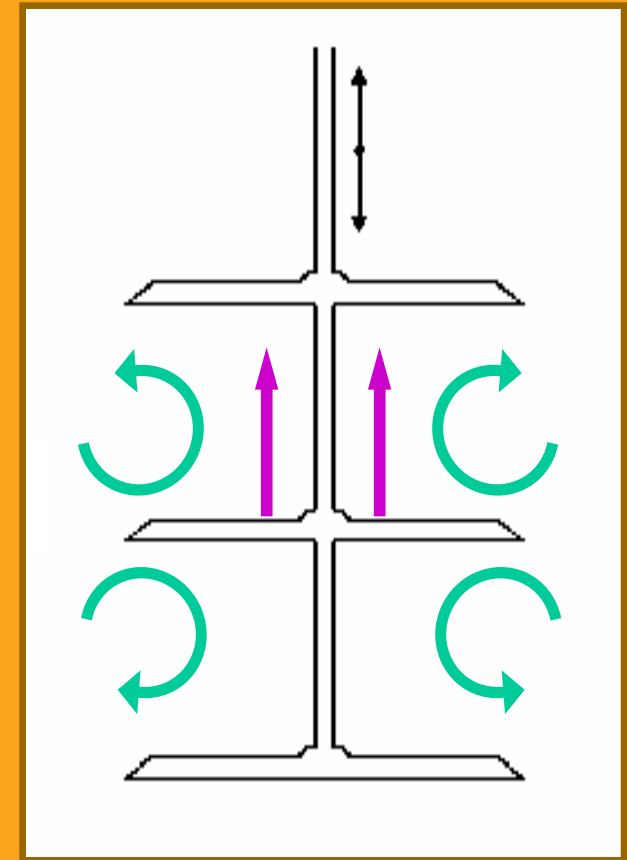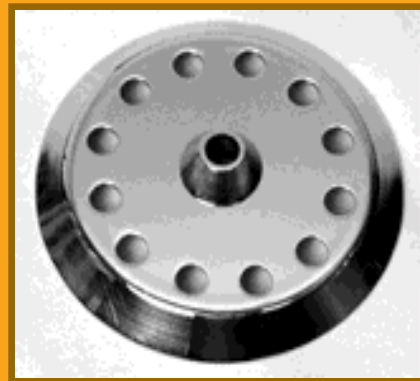

# NEW STIRRING SCHEME

## Stirring properties

- Whole volume stirring
- Good gas distribution
- No vortex formed
- No baffles needed
- Reduced foaming
- Gentle to cells






# NEW STIRRING SCHEME

In action...




# AIR SPARGERS

## a) Capillary

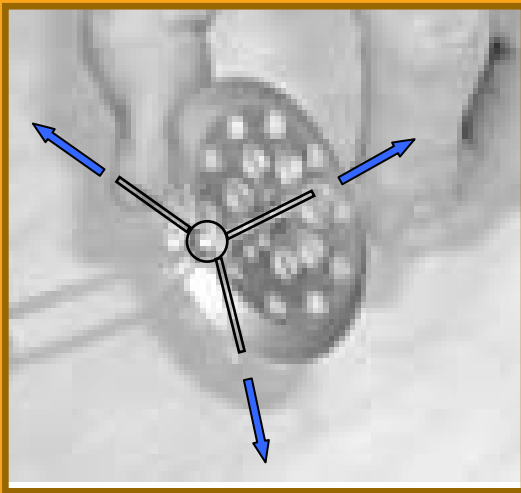

Possible blocking by salts


## b) Silicone

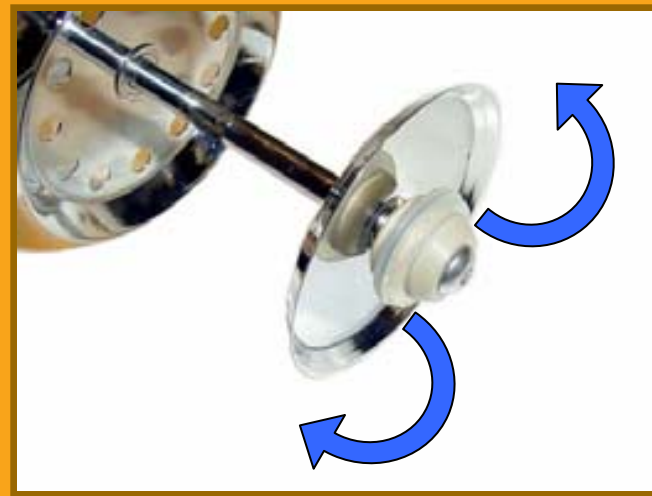

- No blocking by salts
- Good bottom circulation and bubble breaking


# FERMENTATION VESSELS

- Adopted from cell culture
- Up to 8 threaded side necks
- Good accessibility
- Variability
- Low cost
- “jacketable”

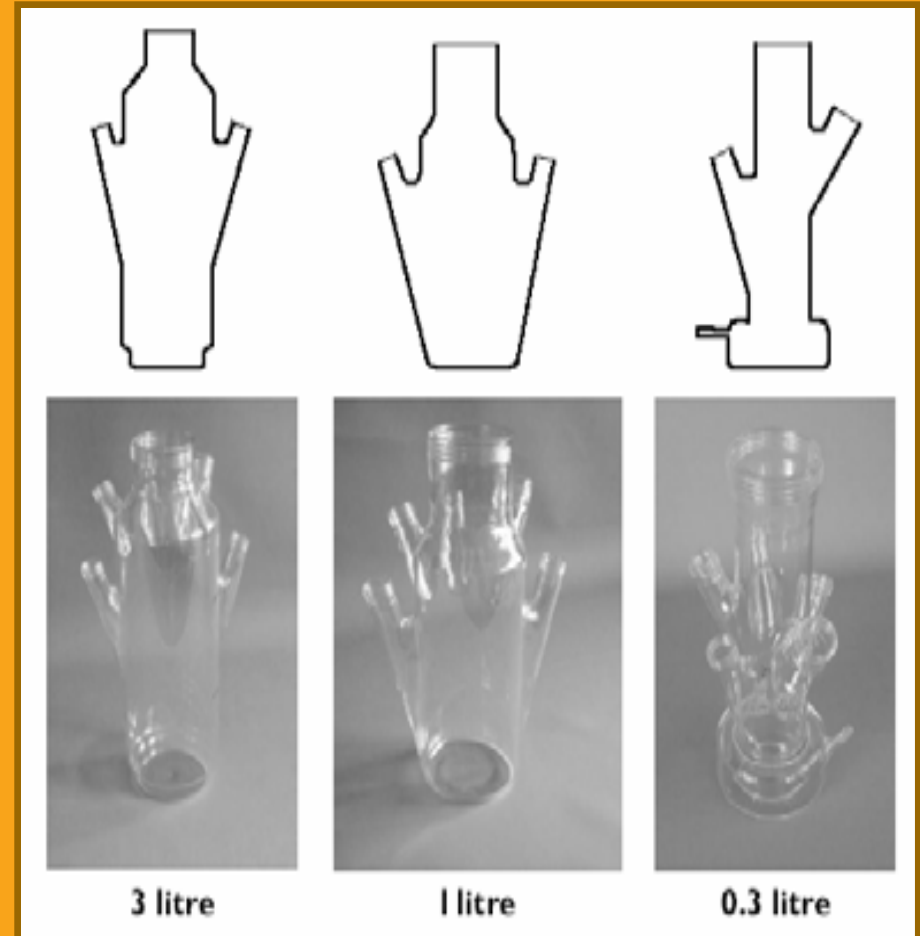

# STOPPER

- Efficient sealing
- Reusable many times

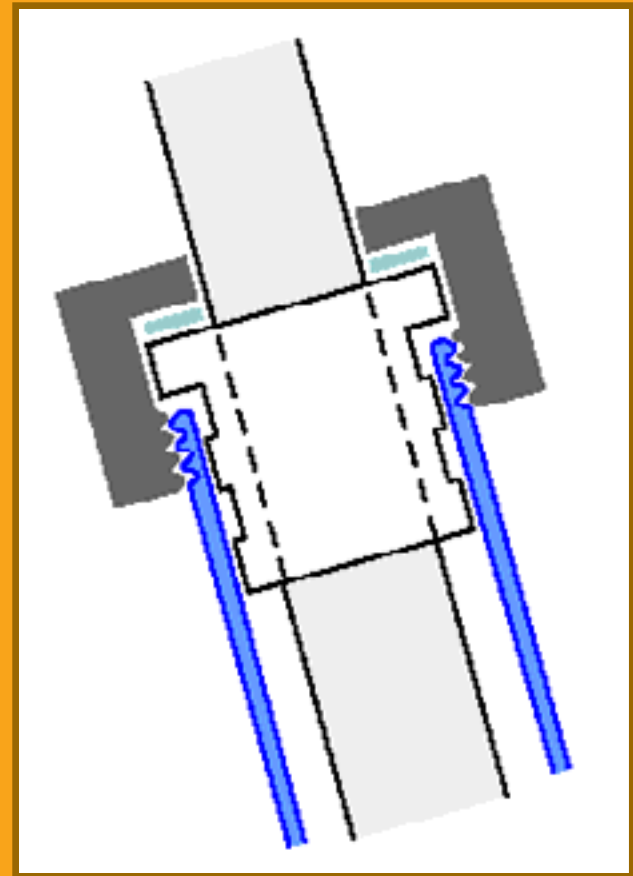

# "COFFEE CUP" RADIATOR

- IR-heating  $\Rightarrow$  gentle sun-like heating
- Natural convection
- No contact with vessel
- No overheating at any volume
- Up to 60°C in 3l vessel
- 98% energy yield
- Negligible heat capacity
- Low voltage



# THE MINIFOR APPROACH

## Fermentor case

- Integrated  $\Leftrightarrow$  free-standing vessel
- Cascaded design
- Ergonomic arrangement
- Excellent accessibility
- Extremely compact



# THE MINIFOR APPROACH

## Control panel

- Closest to operator
- All information at a glance (no scrolling or menu selection)
- Easy operation





# THE MINIFOR APPROACH

## Radiator

- Sealing vessel holder  
⇒ spill protection
- Vessels above heat source
- Shielding of the hot coil  
⇒ protection against burning





# THE MINIFOR APPROACH

## Front connectors

- Shortest connection between probes and vessel
- Sealed cables



# THE MINIFOR APPROACH

## Fermentor platform

- Air output and storage bottles behind the vessel
- Shortest connection between pumps and vessel
- Magnetic holders



# THE MINIFOR APPROACH

Rear side



# THE MINIFOR APPROACH

## General view

- Up to 4 pumps on supports
- Pump protection in case of tubing rupture
- Good visibility of pump status
- Easy accessibility from all sides



# THE MINIFOR APPROACH

## Serial fermentation



- Optimisation in series
- Can be used separately (completely independent)

# PROBES

## pH and temperature probe



- Variopin connector (freely sterilisable)
- Pt 100 integrated in the glass electrode

# PROBES

## DO probe



- PEEK body
- Freely sterilisable connector
- Large cathode system



# PROBES

## DO probe membrane module

- Glass supported PTFE membrane
- Fast response
- Membrane protection





# INOCULATION & SAMPLING

- Stainless steel capillaries
- Luer-Lock adapters



# PERISTALTIC PUMPS



- Constant compression control
- Large rollers
- Asymmetric head
- Compact
- Reliable
- Economy of tubing
- Progressive speed control

# PERISTALTIC PUMPS

In action...



# MASSFLOW

## Gas flow control

- Controlled automatic addition of CO<sub>2</sub> or N<sub>2</sub> during cell cultures
- No need of expensive gas stations
- Programmable
- Totaliser



# CELL CULTURE ADAPTOR




- Very gentle stirring
- Oxygenation through moving coil

# I NTEGRATOR

- Totalisation of delivered liquid or gas
- Access to growth kinetics
- Enzyme activity measurement




# FERMENTATION

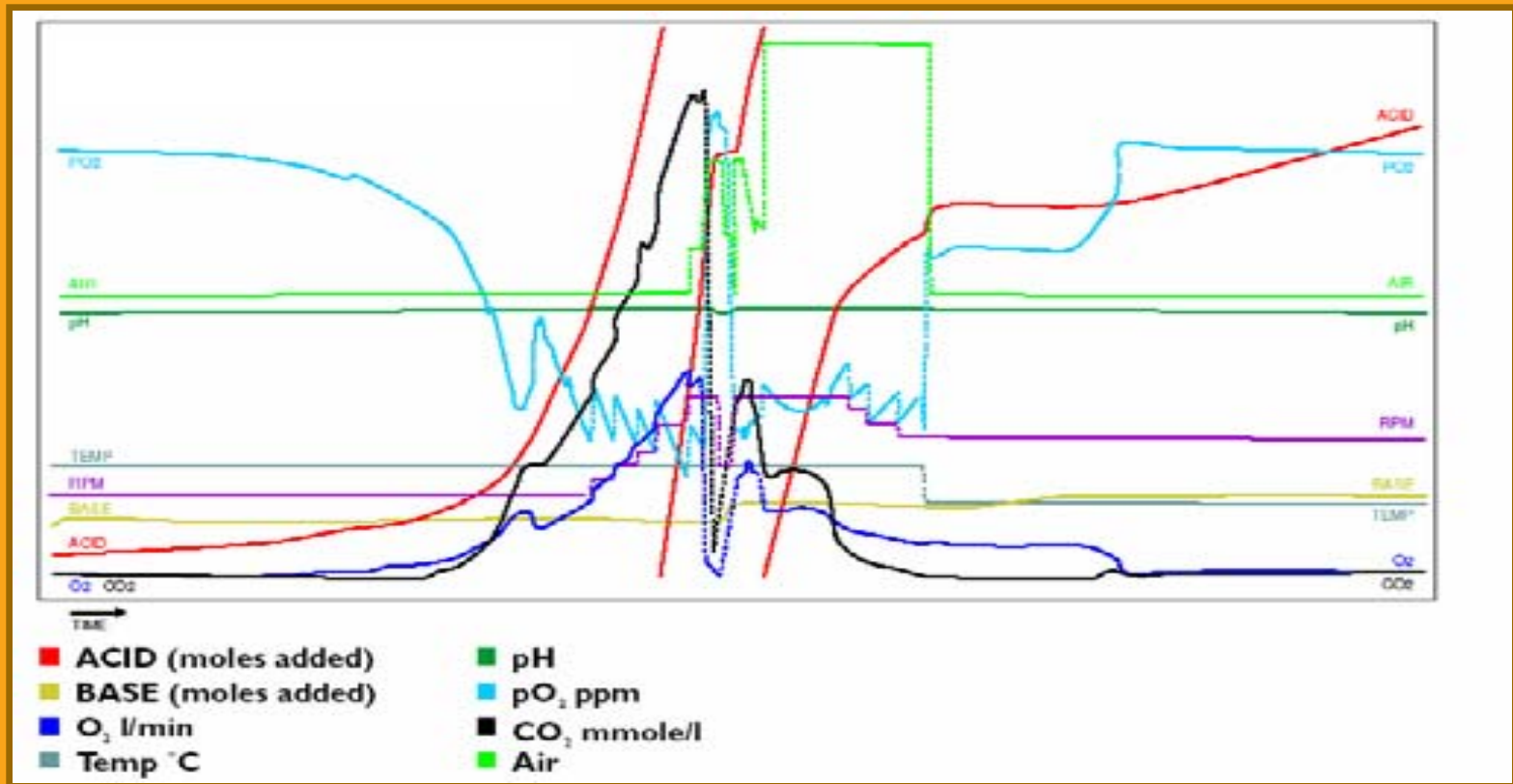

Addition of acid or base necessary to keep the pH constant would not appear without PRECI FLOW and INTEGRATOR !

# FNet - SOFTWARE (Sysmateg)

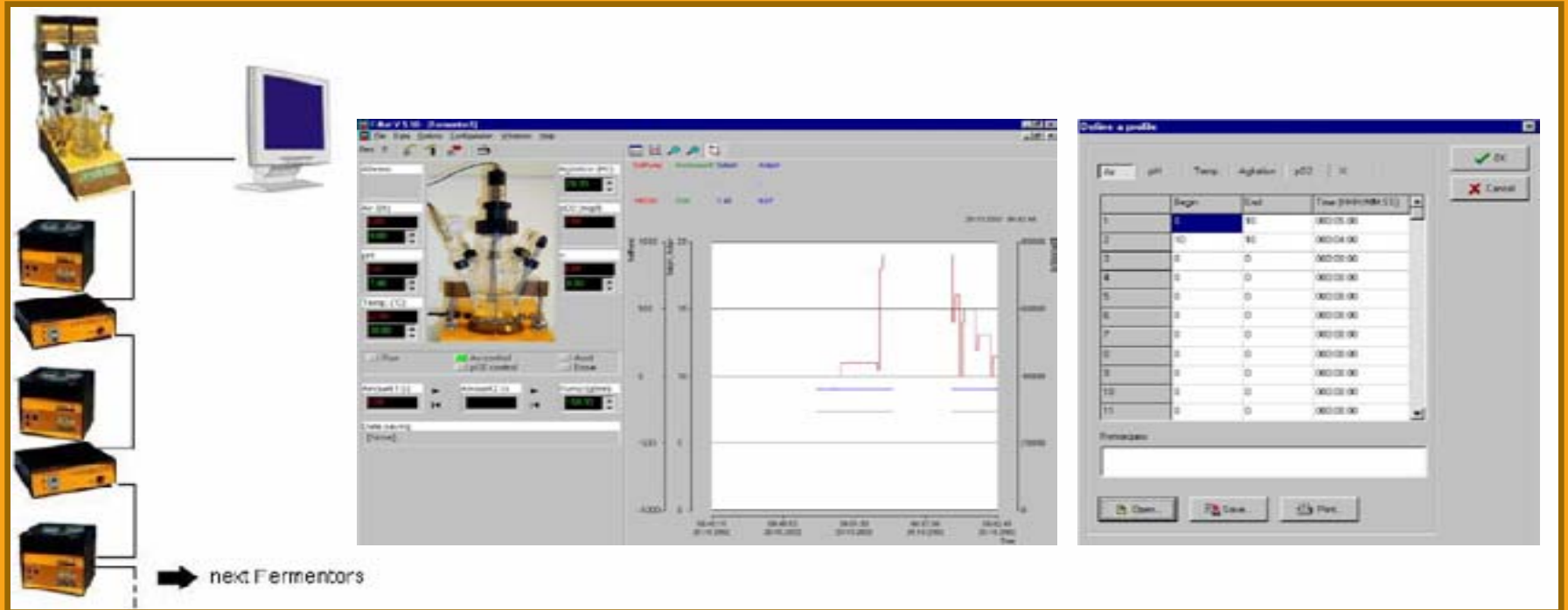

- Complete easy to use software for laboratory fermentors
- Connection of up to 6 fermentors, 12 integrators and 6 pumps

- Data saving
- Graph scaling
- Profile definition
- Alarms



# CONCLUSIONS

- A major effort was invested to make bioreactors more efficient and simple
- New up and down stirring concept providing easy sterility
- IR-heater for gentle heating
- pH and DO probes
- Pumps and gas flow controller

**This work would not have been  
possible without:**

**Jaroslav Bernášek, Vladimír Čermák,  
Rainer Glöckler, Vlastimil Hanuš, Hans  
Kulla, Lida Lehká, Jan Marc Lehky, Petr  
Lehky, Stanislav Tříška, Daniel Venetz,  
Patrick Yax, Michal Zemek**

**THANK YOU !!!**